

MULTI-PURPOSE DECOLORED SHAKER MODEL TS-2000A

MULTI-PURPOSE DECOLORED SHAKER MODEL TS-2000A uses the advanced technology and new shape, It can be with multi-purpose. The upper plate is made from the latex of resisting erosion, the shaker rack can be easily assembled and unassembled, it can mix 2pcs test tube racks which has 4pcs 250ml conical flasks. The accessories can be easily assembled and unassembled, its operation is very simple. The machine is with brushless motor of closed cycle, stepless adjust speed, low noise, without maintain. The machine is with 120 minutes timing and the customer can choose the shaking time and speed.

MULTI-PURPOSE DECOLORED SHAKER MODEL TS-2000A is the equipment for shaking. For decolored: It can be used for the fix of electrophoresis gelatin, the shaking of coloration and decoloration, and It is also used the fix, coloration, developing when the silver nitrate is decolored, and X-ray developing, fixation of autoradiography; and future process of cellulose film electrophoresis; and react and coloration of antigen- antibody. When it is assembled shaker rack, it can be used for the culture for cell and animalcule, and also for the experimentation and study of kinds of shaking, mix and culture.

Main technical data:

Frequency: 40~240 rpm (stepless adjust speed);

Timing: 0~120 minutes;

Power: 85W;

Swing: semidiameter 15mm;

Power supply: AC 200V 50Hz;

Circumstance: -10 degree ~ 45 degree.

Usage:

1. Unpack it, and take the fitting paper under plate, and put the power supply plug into the behind of machine, and put through the power supply, and clean the plate;
2. Put the container of sample on the plate, please note, it should be put placidly;
3. Turn on switch of power supply, the red indicator light is on, and adjust the timing knob to the time you need, and adjust the speed knob from small to big to find the most suitable working.
4. The timing is 0~ 120 minutes, it can be work continuous. If you need continuous work, please adjust the knob into the blank under 0~120, the blank means continuous work.
5. When finish, please turn off the power supply, take away the sample, clean the plate to next operation.

Note: the shaker rack is optional.